MODELL 1 - 10

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MODELS 1 - 10

Our safety cabinets for the storage of hazardous substances are provided with a DN 80 exhaust air connection. The forced ventilation can be carried out by one of our standard extraction fans or by a suitable in-house air extraction system. The exhaust air is to be discharged from the building into a hazard-free zone.

The operation of safety cabinets for the storage of flammable liquids as per DIN EN 14470-1 is governed by TRGS 510. When operating the cabinets without technical forced ventilation, additional replacement measures have to be taken to ensure comparable explosion protection. The herfore required risk assessment we suggest to undertake as per BGR 104 (EX-RL) example collection 2.2.8.

For this reason, and also to avoid the presence of hazardous vapors in the atmosphere near the cabinet, we always recommend that safety cabinets are operated with an extraction fan (min. air exchange rate 10 times per hour) for proper operation in accordance with TRGS 510, annex 1! If an extraction fan is to be used in an EX-endangered zone then it is essential that an EX-proof version is used!

Models 1 + 3 + 9

Extraction from individual safety cabinets

For mounting directly on the storeLAB® safety cabinet - complicated connections from the extraction fan to the safety cabinet are no longer necessary!!

- Not suitable for fume hood underbenches
- With mounting plate for easy mounting on cabinet roof - suction side of fan is connected

If the safety cabinet is used together with an extraction fan then it is no longer necessary to set up and maintain an EX-zone around the safety cabinet!

Installation example: Models 1 + 3 + 9

Models 5 + 7 + 10

Extraction from up to 2 safety cabinets

Extraction fan with mounting base, for separate installation beside (e.g. on wall of room) or on the safety cabinet

- Suitable for all storeLAB® safety cabinets
- With mounting plate for easy mounting on cabinet roof

Installation example: Models 5 + 7 + 10

Models 4 + 8

Plug-in versions

Easy extraction fan connection (suction side) with DN 80 flexible aluminum tubing, with earthed plug and 2 m connection cable.

- Not suitable for fume hood underbenches
- For mounting on safety cabinet

Model 8 with exhaust air monitoring:

- Extraction fan operation checked by monitoring the air flow
- Acoustic message if extraction fan stops working or safety cabinet ventilation slots become blocked

Installation example: Model 8

Description	Model 1	Model 3	Model 9	Model 5	Model 7	Model 10	Model 4	Model 8
EX-proof	NO	YES	YES	NO	YES	YES	NO	NO
Electricity supply	230 V	400 V	230 V	230 V	400 V	230 V	230 V	230 V
Weight (kg)	6,0	6,0	6,0	6,0	6,0	6,0	7,5	7,5
Article no.	B80-2839-A	B80-2841-A	B80-2847-A	B80-2844-A	B80-2846-A	B80-2848-A	B80-2842-A	B80-2843-A

All extraction fans comply with EG Directive 94/9/EG (ATEX)

Connection work to be done by customer!